
The information below is designed to assist you in selecting a powder hood for the ProFiller 3700 or 3800. If you have further questions, feel free to contact CapsuleConnection.com

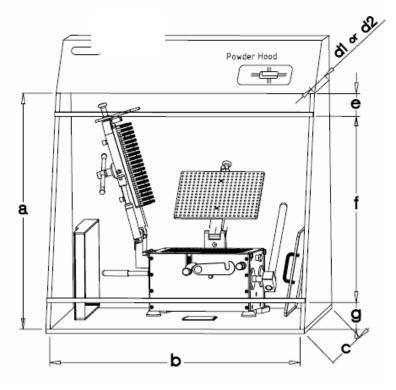

Outside the Powder Hood:

- Empty capsules and filled capsules.
- Orienter.
- Capsule orientation.
- Caps Tray when not in use for separation or locking
- Inside the Powder Hood:
- Filler
- Vibrator (if used)
- Powder Tray¹, Push Plate¹, Powder Spreader
- Caps Tray for separation and locking.

Powder Hood for ProFiller 3700/3800 Filler		Minimum Inner Dimensions Includes Allowance for workspace ²				Top Sash Height ³	Opening Required	Bottom Sash Height ³
		Height	Width	Depth		(optional)	For Operation ³	(optional)
				Bottom	Тор			(
		(a)	(b)	(c)	(d)	(e)	(f)	(g)
Filler	inch	24	27	23	18	6	11 - 15 ⁴	7
	cm	61	68	58	46	15	28 - 38	18

¹ Powder Tray and Push Plate kept vertically against side walls of powder hood.

² The allowance for workspace is 2 inches (5cm) on the height, 2 inches (5cm) on each of the left and right side (width) and 1.5 inches (4cm) on each of the front and back sides (depth).

³ The dimensions "e", "f" and "g" are given considering a table height of 32 inches (73 cm). Dimensions may change based on the height of table. Recommended table height is 29-34 inches (73-86cm).

⁴ Recommended Opening Required for Operation is 11-15 inches (28-38 cm), 11 inches (28 cm) is functional whereas 15 inches (38 cm) is optimal.